

Bottles, water sampling PET

Water test bottles PET, available plain or dosed with Sodium Thiosulphate to inhibit the action of Chlorine, Bromine and Ozone dissolved in water, filled to the requirement of ISO 19458 at 20mg per litre.

All the bottles carry a printed label, have a tamper evident white screw cap and have been sterilised by gamma irradiation at between 25-40 kgy.

Cat. No.	Description	Pk Qty
Undosed		
BO/80000	Bottle PET, water sampling, labelled, sterile, undosed, 330ml	54
BO/80002	Bottle PET, water sampling, labelled, sterile, undosed, 500ml	40
BO/80004	Bottle PET, water sampling, labelled, sterile, undosed, 1000ml	18
Dosed with Sodium Thiosulphate		
BO/80020	Bottle PET, water sampling, labelled, sterile, dosed, 330ml	54
BO/80022	Bottle PET, water sampling, labelled, sterile, dosed, 500ml	40
BO/80024	Bottle PET, water sampling, labelled, sterile, dosed, 1000ml	18



BO/80000

Flocculators, water testing

Flocculator, SW6

For flocculation (jar) tests on water and effluent samples. The SW6 accommodates up to six samples and is specially designed for repeatable conditions between samples and from run to run.

A digital display clearly indicates the speed of the rotational stainless steel paddles, which can be varied from 25 to 250rpm. The digital timer can be set to count down from 1 to 99 minutes.

After count down, the paddles stop and an alert sounds. The timer then counts up in minutes in order to measure settling time. Two pre-set programs allow routine speed and time parameters to be input and then selected by simply pressing one button. Parameters are stored in the memory and can be recalled or altered at any time. The intuitive touch sensitive keypad is totally flat for easy cleaning. Sample monitoring is assisted by diffused base illumination and black background.

A simultaneous coagulant doser is available as an accessory.

- Accepts six samples
- Digital indication of speed and time
- Two pre-set programs
- Built in illumination for easy viewing of samples

Flocculator, SW5

A mini version of the SW6 that accepts two samples (without timer function). The case incorporates a handle for easy transportation.



WT/15002

Technical Specifications		
	SW5	SW6
No. of samples	2	6
Speed range	25 to 250rpm	25 to 250rpm
Timer	No	0 to 99 minutes
Pre-set programs	2	2
Controls keypad	Dial	Touch sensitive
Digital displays	LCD	LED
Dimensions (WxDxH)	320 x 210 x 500mm	750 x 210 x 460mm
Net weight	8.6kg	17kg
Power	230V, 50Hz, 100W	230V, 50Hz, 100W

Cat. No.	Code	Description
WT/15002	SW6	Flocculator with 6 rotators (without beakers)
WT/15004	SW5	Flocculator with 2 rotators (without beakers)
WT/15006	SW6/2	Simultaneous coagulant doser (SW6 only)
BE/12276	1015/22D	Pyrex® glass beakers, 1000ml, pack of 10



WT/15004